

# WGLT

## Wireless Glove Leak Tester

### Introduction

The **Esco Wireless Glove Leak Tester (WGLT)** facilitates the detection of any punctures or holes to guarantee the integrity of an isolator's gloves/sleeves system. This accessory uses the principle of pressure loss and method based on Class 3 Containment as per ISO 10648-2.

### The WGLT comprises of two parts:

- 1. Hardware** - refers to the wireless glove leak tester and its independent control.
- 2. Software** - refers to the application used to remotely control and monitor the glove leak tester. This is generally installed on a Windows-based computer which meets the system requirements.



### Quick View



1. Seal Deflate Button

2. Control Button with Operation Status LED

3. USB Port

4. Power Button with Battery Status LED

5. Charging Port

For items #2 and #4: The button will light up to showcase the WGLT's running status. Refer to the color legend below the buttons.

### Key Features





- Sleek, easy to clean, and low-maintenance design
- FDA-approved inflatable seals
- HDPE glove port disc for increased durability
- Portable and wireless for on-the-go use
- Generates alarm and data logs and can be operated within assigned authority levels
- Remotely connects and controls the unit
- Glove port docking part can be customizable to match any existing glove port size. (Note: Actual sample of the glove port will be needed for accurate sizing of the WGLT)

### Basic Principles

- Inflatable seals inflate to securely seal the glove
- The compact built-in pump of the glove leak tester injects air into the glove until it reaches the desired pressure
- The glove leak tester then stops the air injection and reads pressure loss within 5 minutes
- The glove leak tester determines whether the glove passes the leak tightness test or not in order to monitor the integrity of the isolator's gloves/sleeves system

SOFTWARE GENERAL SPECIFICATIONS	
Component	Recommended
Processor (CPU)	Intel Core i5 or equivalent
Operating System	Windows 10 Professional x64
Memory (RAM)	8 GB
Network Adapter	802.11AC Dual Band 2.4g/5g GHz wireless network adapter
Others	.NET Framework 4.6.1, MS Access Database Engine, Crystal Report Runtime (included in the installer package) Router that supports IPv4 NAT or IPv6 Drive storage consumption: 34.8 MB











HARDWARE GENERAL SPECIFICATIONS	
Glove type	Oval Glove
External Dimension	295 mm x 189 mm x 395 mm
Net Weight	3.75 kg
Disc size	208 mm x 308 mm
Exterior Material	HDPE
Battery	14.8 V, 5 AH (21700 Li-Ion Battery 4S1P)
Charger	16.8 VDC, 1.8 A; 3 hours computed charging time
Number of Tests	8 full cycle tests
Maximum pump pressure	250 kPa
Pressure Range	0 – 300 Pa
Processor	Single Core ARM 1 GHz
RAM, Runtime memory	512 MB
ROM, Data, and Log Memory	8 GB Flash MicroSD
USB Interface	1 x USB 2.0 Type A

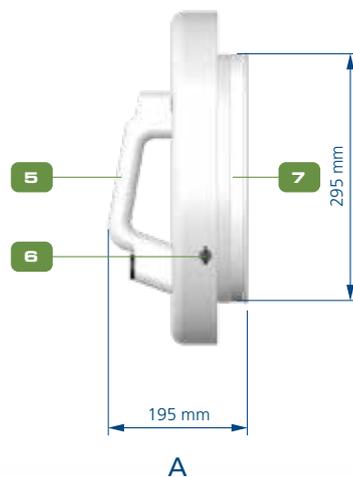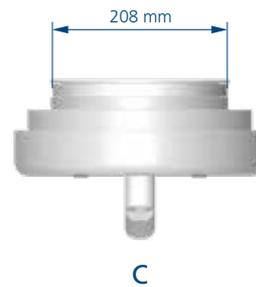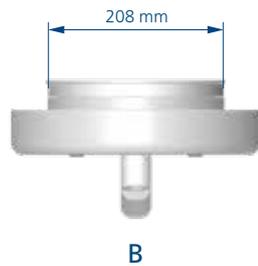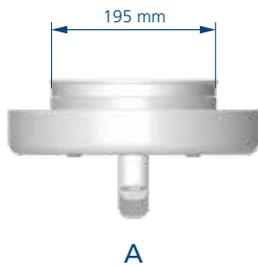

1. Seal Deflate Button
2. Control Button with Operation Status LED

3. USB Port
4. Power Button with Battery Status LED

5. WGLT Handle
6. Charging Port
7. Inflatable Seal

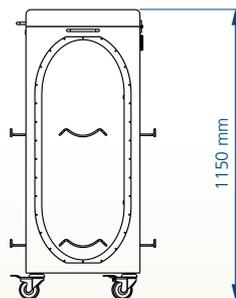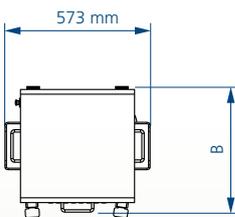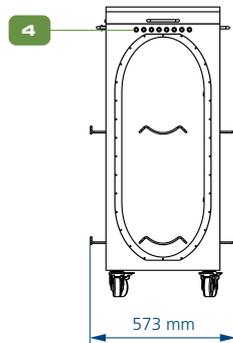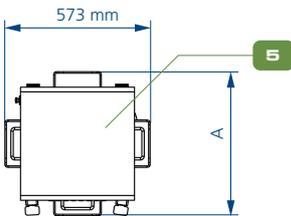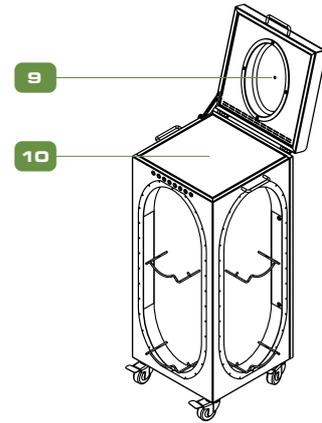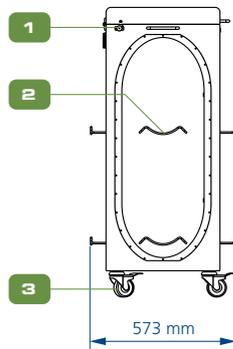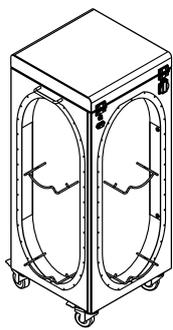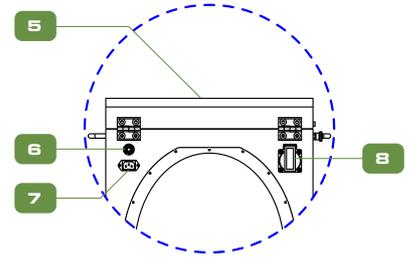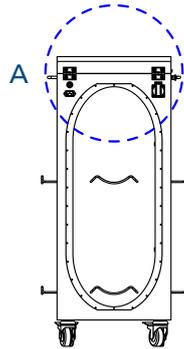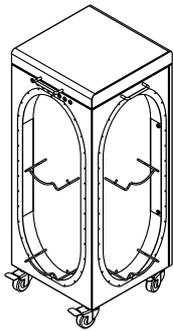

1. Retractable Spring Plunger
2. WGLT-G2 Holder
3. Castor Wheel With Brake
4. Push Button Switches For WGLT-G2 Charger
5. Hinged Outer Table
6. Circuit Breaker Reset Button
7. Power Supply Inlet
8. Electrical Socket IP54, Type G
9. Inner Table
10. Test Port

# ESCO LIFESCIENCES GROUP NETWORK

42 Locations in 21 Countries All Over the World



- Air Shower
- Aseptic Containment Isolator (ACTI)
- Ceiling Laminar Airflow Units
- Cleanroom Transfer Hatch
- Containment Barrier Isolator (CBI)
- Downflow Booth (DFB)
- Dynamic Floor Laminar Hatch
- Dynamic Pass Box
- Evidence Drying Cabinet
- Garment Storage Cabinet
- General Processing Platform Isolator (GPPI)
- Laminar Flow Horizontal Trolley
- Laminar Flow Straddle Units, Single and Double
- Laminar Flow Vertical Trolley
- Pass Box
- Soft Wall Cleanroom
- Sputum Booth
- Ventilated Balance Enclosure (VBE)
- Weighing and Dispensing Containment Isolator (WDCI)

Since 1978, Esco has emerged as a leader in the development of controlled environment, laboratory and pharmaceutical equipment solutions. Products sold in more than 100 countries include biological safety cabinets, fume hoods, ductless fume hoods, laminar flow clean benches, animal containment workstations, cytotoxic cabinets, hospital pharmacy isolators, and PCR cabinets and instrumentation. With the most extensive product line in the industry, Esco has passed more tests, in more languages, for more certifications, throughout more countries than any biosafety cabinet manufacturer in the world. Esco remains dedicated to delivering innovative solutions for the clinical, life science, research and industrial laboratory community. [www.escoglobal.com](http://www.escoglobal.com).

### Esco Pharma

21 Changi South Street 1 Singapore 486777  
 Tel: +65 65420833  
 Email: [csis.pharma@escolifesciences.com](mailto:csis.pharma@escolifesciences.com)

### Esco Technologies, Inc.

2512 Metropolitan Drive, Suite 120 B  
 Feasterville- Trevose, PA 19053-6738  
 Tel: +1 215 322 2155  
 Email: [eti.pharma@escolifesciences.com](mailto:eti.pharma@escolifesciences.com)

### Esco GB Ltd

Unit 2 R-evolution @ Gateway 36, Kestrel Way, Barnsley, S70 5SZ  
 Tel: +44 (0) 1226 360 799 • Email: [egb.info@escolifesciences.com](mailto:egb.info@escolifesciences.com)

**Esco Lifesciences Offices:** Bangladesh | China | Denmark | Germany | Hong Kong | Indonesia | Italy | Lithuania | Malaysia | Myanmar | Philippines | Russia | Singapore | South Africa | South Korea | Taiwan | Thailand | UAE | UK | USA | Vietnam



Wireless\_GLT\_sellsheet\_A4\_vC\_07102023  
 Esco can accept no responsibility for possible errors in catalogues, brochures and other printed materials. Esco reserves the right to alter its products and specifications without notice. All trademarks and logos in this material are the property of Esco and the respective companies.

